


Resetting the over temperature cut-out

The over temperature cut-out on the heater mat tripping out is generally caused by the water bath trying to heat with no or very little water in it or by large objects being placed directly on the bottom of the tank with no tray in place.

To reset the over temperature cut-out first remove all power and the water from the water bath. Then turn the unit upside down and remove the bottom cover, by removing the four or six screws (depending on the size of bath).

Note: the feet are fitted to the bottom cover and do not require removing.

The bottom of the tank is now visible and on that is the heater mat (see below), in the middle of the heater mat is the cut-out. In the centre of the cut-out is a button this requires to be pressed in.

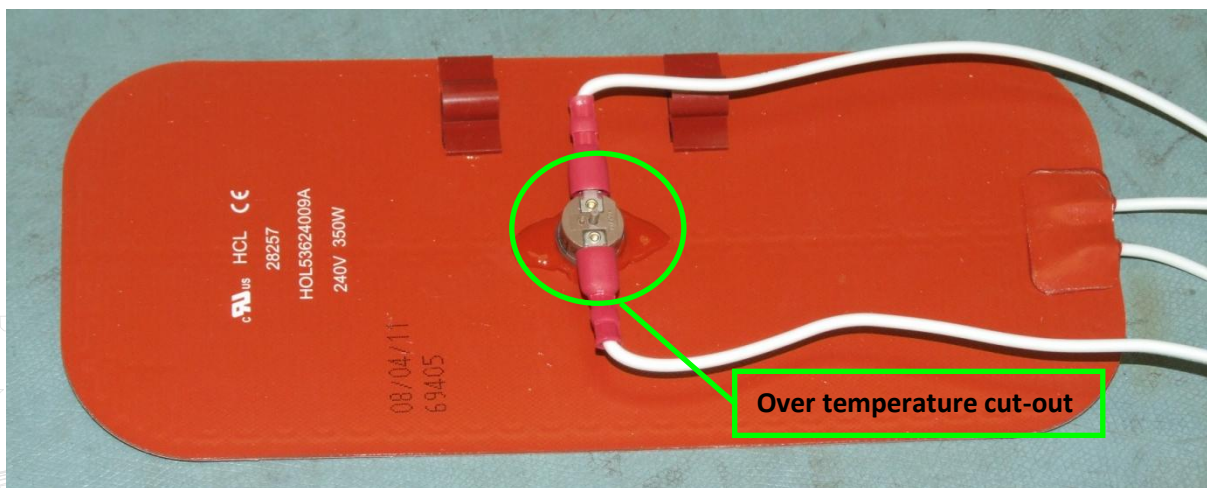

Refit the bottom cover and fit the original four or six screws removed.

If the cut-out has been tripped this will reset the bath and restore to normal operation. Ensure that the trip cause is known and corrected before further operation.

Reconnect the water bath to the power and refill with water before turning it on.

If you require further information please contact your local distributor or Grant Instruments technical support department at scientifictech@grantinstruments.com